

VS Suspension Solution

rev.4/25

CAT#:	TCR-SS001-100
AMOUNT:	100 mL
FORM:	Liquid
STORAGE CONDITIONS:	4°C
SHIPPING:	Room Temperature
STERILITY:	0.22 µm Sterile filtered under Sterile conditions
SHELF LIFE:	Stable for 12 months when stored and handled under proper conditions

DESCRIPTION:

VS Suspension Solution is a protein free buffer solution optimized for diluting VisionSort calibration and performance beads.

INSTRUCTIONS FOR USE:

- For VS C-Beads, dilute **10µL** of VS C-Beads with **500µL** of VS Suspension Solution.
- For VS P-RBW Beads, dilute **20 µL** of VS P-RBW Beads with **500 µL** of VS Suspension Solution.
- For Vision Beads, dilute **120 µL** of Vision Beads with **120 µL** of VS Suspension Solution.

UNITED STATES

1100 Island Drive, Suite 203, Redwood City, CA

JAPAN

7-3-1 Hongo, Bunkyo, Tokyo

THINKCYTE.COM

Copyright © 2024 ThinkCyte, K.K.. ThinkCyte, Ghost Cytometry, and VisionSort are trademarks of ThinkCyte, K.K.. All rights reserved.
For Research Use Only.